

[www.cranepharmaceutical.com](http://www.cranepharmaceutical.com)

Saunders<sup>®</sup> Performance Manual Bonnet and EC Compact actuator are both manufactured from PES (Polyethersulphone), a high performance thermoplastic. PES has outstanding strength, chemical resistance and thermal properties to maximize product integrity.

The chemical resistance of PES has been investigated and a summary of the resistance of PES to various chemicals is provided (on the back page). The results are derived from tests in which unstressed specimens were completely immersed in a wide range of chemical environments at room temperature.

In common with most thermoplastics, the resistance of PES to chemicals is also dependent on stress levels and on temperature. These conditions, particularly those of stress and strain, are difficult to reproduce in the laboratory. The table, therefore, should be used as a guide, and the user should satisfy themselves beforehand of the suitability of PES for the in-service environment.



PES is also used as standard hand wheel material for Saunders<sup>®</sup> Stainless Steel Rising Hand Wheel Bonnet and Performance Bonnet.

## PES Chemical Resistance Summary

A summary of the resistances of PES to various chemicals is given below:

### Water

PES is not attacked chemically by water.

### Inorganic Chemicals

PES is not unaffected by most inorganic reagents. It has good resistance to aqueous acids, bases, and most inorganic solutions. PES is attacked by concentrated oxidising mineral acids (e.g. Nitric and Sulphonic Acids) at room temperature but is not affected by more dilute acids. Resistance to Alkalis is good.

### Sterilizing Solutions and Anaesthetics

Most of the common sterilizing solutions and anaesthetics can be used safely in contact with PES at their recommended concentrations.

The excellent resistance of PES to hot water means that it can be repeatedly steam sterilized. PES can also be sterilized by radiation as it offers high resistance to beta gamma and x-rays.

### Halogenated Cleaning Solvents

Many cleaning and degreasing solvents are based on chlorinated and fluorinated hydrocarbons. Unless heavily stressed, PES can be cleaned by most of these solvents.

Solvents not recommended are methylene chloride, 1,1,2-trichloroethane or chloroform. (Some proprietary cleaning solvents consist of 1,1,2-trichloroethane diluted with other solvents - these are suitable for PES.)

### Aromatic Solvents

Components made from PES can be used for short-term exposure to solvents such as benzene, xylene, and toluene, provided they are not heavily stressed. Polar aromatic solvents, such as N-Methyl Pyrrolidone, are often solvents for PES. Overall avoid contact with these types of solvents.

### Oil and Greases

Test show that PES is suitable for prolonged contact with most oils and greases at high temperature. The only known exceptions are di-ester and phosphate-ester based oils.

### Petrol

At room temperature, PES can be used in contact with Petrol.

### Solubility Properties of PES

PES is a polar polymer and therefore tends to dissolve in polar solvents, for example:

- Dimethyl Sulphoxide
- N, N-Dimethylformamide
- N-Methylpyrrolidone
- Pyridine
- Quinoline
- Aniline
- O-Chlorophenol

## The resistance of PES to Chemical Attack

A = No attack. Little or no absorption at 20°C (68°F)

B = Slight attack. Some absorption causes swelling at 20°C (68°F.) Satisfactory use of PES will depend on the application.

C = Bad attack. PES should not be used for any application where these environmental conditions are present.

The use of a small letter indicates that PES has not been tested specifically with that chemical but that the result can be predicted from tests with similar chemicals.

Chemical	Result	Chemical	Result	Chemical	Result
Acetaldehyde	C	Dichlorobenzene	C	Nitrobenzene	C
Acetic Acid - Glacial	A	Diesel Oil	A	Oils - Vegetable	A
Acetic Acid - 10%	A	Diethylamine	A	Oleic Acid	a
Acetone	C	Diethyl Ether	A	Oleum	C
Aluminium Salts	a	Dimethyl Formamide	C	Oxalic Acid	A
Ammonia 880	A	Diocetyl Phthalate	A	Perchloroethylene	C
Ammonium Hydroxide - 10%	A	Dioxane	b	Petrol	A
Ammonium Chloride - 10%	A	Edible Fats and Oils	A	Petroleum Ether	A
Amyl Acetate	B	Ethyl Acetate	C	Phenol	C
Aniline	C	Ethyl Alcohol	A	Potassium Hydroxide - 10%	A
Arcton Propellants	a	Ethylene Glycol	A	Potassium Hydroxide - 50%	A
Aviation Hydraulic Fuels	B	Ferric Chloride	A	Propane	a
Aviation Spirit	A	Formaldehyde	A	Pyridine	C
Barium Salts	a	Formic Acid	A	Silicone Fluids	A
Benzaldehyde	C	Glycerol	A	Silver Nitrate	A
Benzene	A	Heptane	A	Soap Solution	A
Benzoic Acid	A	Hexane	A	Sodium Chloride	A
Benzene Sulphonic Acid	a	Hydrochloric Acid 10%	A	Sodium Hydroxide - 10%	A
Bleach	A	Hydrochloric Acid Concentrate	A	Sodium Hydroxide - 50%	A
Boric Acid	a	Hydrogen Peroxide	A	Sodium Hypochlorite	a
Brake Fluid	B	Hydrogen Sulphide	a	Sulphuric Acid - 10%	A
Brine	A	Iodine in Potassium Iodide	B	Sulphuric Acid - Concentrate	C
Butane	A	Isopropanol	A	Sulphurous Acid	c
Butanol	A	Iso-Octane	A	Tar	a
Butyl Acetate	b	Kerosene	A	Tartaric Acid	a
Calcium Nitrate	A	Lactic Acid	a	Tetrahydrofuran	C
Calcium Hypochlorite	A	Lead Acetate	a	Toluene	C
Carbon Disulphide	b	Linseed Oil	A	Transformer Oil	A
Carbon Tetrachloride	A	Magnesium Sulphate	A	Trichloroethylene	B
Chlorobenzene	C	Mercuric Chloride	a	Turpentine	A
Chloroform	C	Mercurous Chloride	a	Vaseline	a
Chlorosulphonic Acid	c	Mercury	a	Varnish	A
Chromic Acid	A	Methanol	A	Water	A
Citric Acid	A	Methyl Ethyl Ketone	C	Wax	a
Copper Sulphate	A	Methylene Chloride	C	White Spirit	A
Creosote	A	Milk	A	Wines and Spirits	a
Cresols	C	Motor Oil	A	Xylene	B
Cyclohexane	A	Nickel Salts	a	Zinc Salts	a
Dichloroethane	C	Nitric Acid - 10%	A		
Dichloroethylene	B	Nitric Acid - Concentrate	C		

Please visit our Web-Based Drawing Library at: [www.saundersdrawings.com](http://www.saundersdrawings.com) for current database of drawings in PDF, 2D DWG, and 3D STP formats.

[www.cranepharmaceutical.com](http://www.cranepharmaceutical.com)



CRANE ChemPharma Flow Solutions, Saunders® • 9860 Johnson Road  
Montgomery, Texas USA 77316 • Tel.: +1 936 588 8360 • Fax: +1 936 588 8302

IMPORTANT: CRANE Co. and its subsidiaries are not responsible for the accuracy, compliance, and legality of material contained in this datasheet offered in print, on the company websites, or via any external links, or third party sites.